2D NMR Spectrum Processing with Mnova

Hongjun Zhou Dept of Chem & Biochem UC Santa Barbara

April 09, 2012

General NMR Data Processing

NMR data processing is typically done in the following order:

- Read in raw FID
- Select and apply apodization or window function
- Zero fill (**ZF**) the FID, usually at least doubling the original data size
- Fourier transform (FT)
- Phase spectrum
- Baseline correction

Some of these are done automatically by Mnova when data are read in.

NOTE:

- Phasing means, in most cases, to adjust the spectrum so that peaks across the spectrum are in-phase (absorptive mode) where all peaks point up (or down) with the peak bases look even on both sides.
- Zero-order phase (PH0) is a constant phase across the spectrum and is applied first. Next, set a pivot point for 1st-order phase (PH1) adjustment. The **pivot** should be ideally set to a strong peak on one end of the spectrum. Apply 1storder phase which linearly changes across the spectrum from the "pivot" point.
- Baseline correction should be applied AFTER phasing. If a small phase adjustment is still needed after baseline correction, adjust the phases after baseline correction.

2D NMR Spectrum Processing with Mnova

The two dimensions in a 2D spectrum is specified as t2 or f2 (horizontal) and t1 or f1 (vertical) dimensions, respectively. t1 and t2 refer to the time-domain data; f1 and f2 refer to the frequency domain data after FT of t1 and t2 dimensions. The processing command below is often applied to f2 and f1 (or t2 and t1) separately. Select either **f2** or **f1** from the top menu buttons before applying the processing command.

| 近 式 | 十 | 弐 ・ || ff1 f2 | ℡ ・ | 🎠 斧 ・ 🌾 ・ 🍂 ・ 楡 🤉

2D spectrum should be processed in the following order:

- 1. Open *.fid data
- 2. Select appropriate window functions for t2 and t1 under *Processing->Apodization*
- 3. Decide whether the spectra need phase adjustment
 - 1. For all absolute-value (or magnitude) mode spectra, such as typical COSY and HMBC, **DO NOT** apply phase correction.
 - For phase-sensitive spectra, such as typical HSQC/HMQC/NOESY etc., phase both f2 and f1 dimensions. Apply both zero-order (ph0) and 1st-order (ph1) phases if necessary.
- 4. Apply baseline correction to **f2** and **f1** dimensions.
- Play with other optional processing tricks such as t1 noise reduction or symmetrization, etc. Be cautious that these routines may introduce artifacts or reduce signal/noise. Always compare results with those from simple processing.

When 1st-Order Phase Is Needed

Two spectra after zero-order phase correction



1st-order phase is NOT needed

1st-order phase is needed. Note the linear change in phase from left to right.

2D Processing Adjustment in Mnova



Magnitude-mode COSY Processing

NOTE: No phase correction should be applied.



gCOSY: Initial look after automatic processing

• Vertical stripes are called t1 noise



Check and set *Processing→apodization* function to Sine Square 0.0 along t2 and t1



Apply baseline correction along f1 (vertical)



Raise the intensity threshold with mouse scroll



Turn on **Processing** \rightarrow **Reduce t1 noise.** In this case, the improvement is significant, but be cautious with artifacts. To turn off t1 noise correction, toggle off **Processing** \rightarrow **Reduce t1 noise**



Use *View→Crosshair* (or press *C*) to check symmetry of COSY crosspeaks Hold the left mouse button and drag to show a box



Select *View→2D Plotting Method→Contour Plot* or right mouse button click. Contour drawing on the fly is slower. Do this after data processing.



NOESY Processing: Initial look

- Reference is wrong
- Apodization/window function need to change



Referencing



Select f1 or f2 from menu buttons and change *Processing*→*Apodization* to: Sine Square 90.0 along t2 Sine Bell 90.0 along t1



The 2D peaks will look smoother and the baseline noise is less with the new window function



- Select manual phase correction
- Select f2. Drag pivot line (vertical blue line) to line up with a strong peak on left
- Adjust PHO so that the selected peak is perfectly in phase
- Adjust PH1 so that the peaks on right also come into phase



After PHO adjustment along f2



After PH1 adjustment along f2



- For f1, select a peak near the bottom (or top) as the pivot
- Spectrum after PHO adjustment
- PH1 seems OK. NO linear phase correction necessary.



- Lower threshold so that baseline defects are seen
- Apply Polynomial baseline correction order 2 or 3 to f2 and f1 separately



After f2 baseline correction



After both f2 and f1 baseline correction



Additional processing

- Processing→Symmetrize→COSY-like
- Spectrum quality improves.

NOTE:

 Be careful with the "reduce t1 noise" or symmetrization procedure. Artificial crosspeaks may be created in the process if the background noise is strong.



- For small molecules, NOEs mostly have opposite signs with the diagonal peaks.
- Typical artifacts are:
 - dispersive peaks (with mixed up/down signals) that come from through-bond COSY-correlations. These are suppressed but not completely.
 - exchange crosspeaks having same sign with diagonal peaks



These through-bond COSY-type peaks tend to cancel some or all through-space NOEs because they have opposite signs in some overlapping areas. Here, more COSY-type peaks are seen between the aromatic ring protons.



Exchange crosspeaks in NOESY are shown as in-phase peaks having the same sign as the diagonal peaks but opposite to NOESY crosspeaks. They can result from:

- chemical exchange, such as exchange of ¹H in H₂O with an amide ¹H.
- exchange among multiple conformers where a ¹H has a distinct chemical shift in each conformer.

